


Core Science Facility Vertical Fume Hood Guidelines



Operating Sash Height

Position sash at 9". 18" should be used minimally as the alternate height.

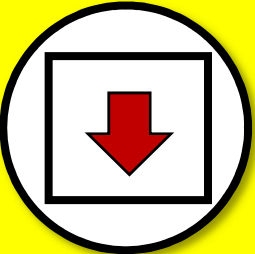

Close Sash

Close sash when fume hood is not actively being used.



Working in Hood

Conduct work at least 6" from sash opening.



Baffle

Ensure baffle is clear of any objects that can obstruct airflow.



Using Equipment

Elevate equipment at least 2" to ensure adequate airflow.



If Alarm Sounds

Verify that sash is closed or is positioned at 9" or 18"
If alarm continues contact techsvs@mun.ca

July 2024



Never disarm or remove fume hood alarm